





# Cellvento® and EX-CELL® Feed Mixing Protocol

1.

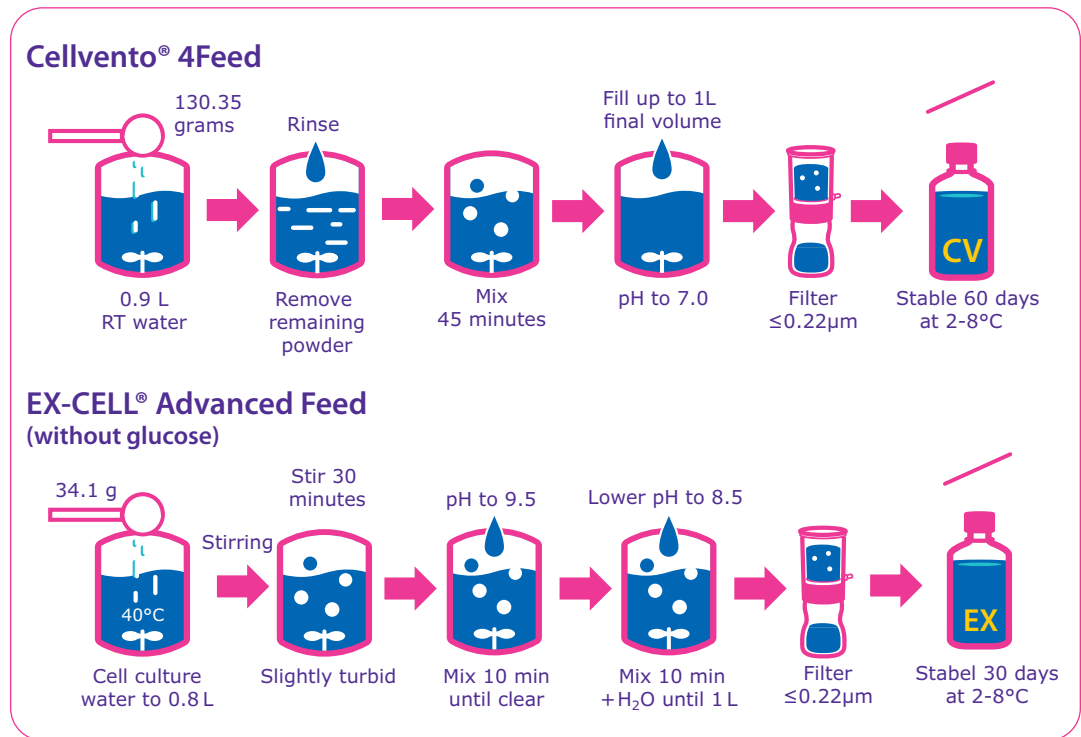

2.

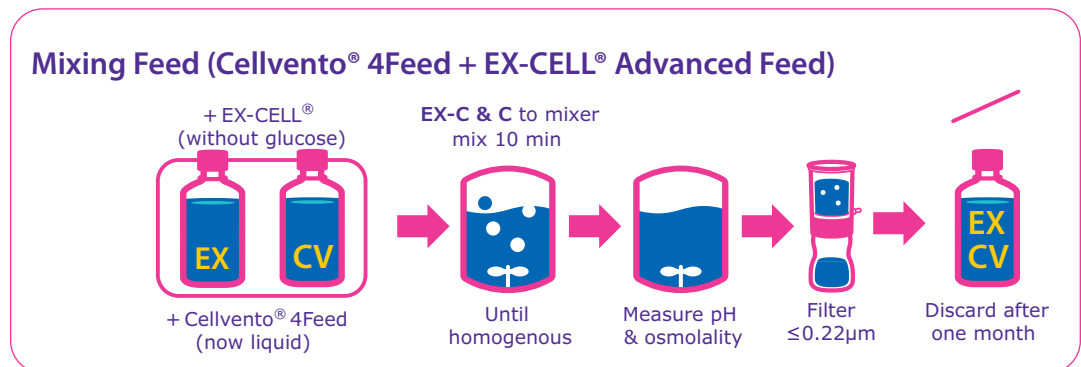

Additional information on back side.

## Try different mix levels to find the optimal level of feed for your process

### Suggested Mixing Levels

Mix Ratio	Cellvento® 4Feed	EX-CELL® Advanced Feed
Volume for one liter of mixed media		
100	1 L	0 L
75:25	750 mL	250 mL
50:50	500 mL	500 mL
25:75	250 mL	750 mL

Evaluate Cellvento® 4Feed as a mix and a standalone feed. Usually EX-CELL® Advanced Feed performs best with the Cellvento® 4Feed.

In some applications, Cellvento® 4Feed may perform better without EX-CELL® Advanced Feed.

### To place an order or receive technical assistance

In the U.S. and Canada, call toll-free  
1-800-645-5476

For other countries across Europe  
and the world, please visit:  
[EMDMillipore.com/offices](https://www.emdmillipore.com/offices)

For Technical Service, please visit: [EMDMillipore.com/techservice](https://www.emdmillipore.com/techservice)

[EMDMillipore.com](https://www.emdmillipore.com)

MilliporeSigma  
400 Summit Drive  
Burlington, MA 01803

