

CellPrime® rAlbumin

Recombinant Albumin

Product Description

CellPrime[®] rAlbumin is a non-animal origin (NAO) recombinant media supplement that is GMP manufactured according to EXCiPACT[™]. It combines excellent quality with superior performance and lot-tolot consistency. This supplement is expressed from a synthetic DNA insert that encodes the human albumin gene sequence in the yeast Pichia pastoris. Like all members of the recombinant CellPrime[®] portfolio, CellPrime[®] rAlbumin is a non-animal origin (NAO) product that does not contain components of animal or human origin in the

- Master Cell Bank (MCB),
- Working Cell Bank (WCB),
- Production raw materials,
- Manufacturing process, or
- Final product

Product Applications

Cellprime[®] rAlbumin does not use octanoate (caprylic acid), a commonly used stabilizer for albumins during pasteurization, in the manufacturing process or in the final product formulation. Octanoate is not well tolerated by a number of cell types.

CellPrime[®] rAlbumin can replace serum albumin (bovine or human serum origin albumin) in combination with recombinant transferrin and recombinant insulin as a supplement mix in cell culture media formulations. It removes the variability introduced into the manufacturing process by blood-derived proteins and safeguards against adventitious virus or prion contaminations. It mitigates the risk of contaminating a production facility that may result in a manufacturing shutdown and hence create supply shortages for patients.



This product aligns with our total viral safety strategy: **Prevent**, Detect, Remove and affords our customers a safe alternative to human or bovine serum derived albumin.

Thus, CellPrime[®] rAlbumin is the preferred supplement for the extension of these cell lines used in either (virusborne) vaccine manufacturing or cell and gene therapies:

- MDCK, Vero, MRC-5, HEK 293 (T)
- Mononuclear and hematopoietic stem cells, T-lymphocytes
- Human mesenchymal stem cells (MSC), human embryonic (hESC) and induced pluripotent stem cells (iPSC)



Key Benefits of Product Offering

>=98% pure, multi-functional protein

- Oncotic pressure regulator
- Carrier for hormones, metabolites, fatty acids, minerals, scavenger of toxic radicals
- Cell therapy and viral vector stabilizer used for formulation and cryopreservation
- Stem and immunotherapy cell culture supplement
- Surface coating agent
- Recombinant, produced without the use of any animal- or human-derived material at any point of the manufacturing process (NAO)
- Recombinant replica of human serum albumin (non-glycosylated)
- Earmarked to attain Emprove® status
- Shelf-life: 2 years at recommended storage temperature

Handling Instructions

- Store and ship CellPrime[®] rAlbumin refrigerated at 2–8 °C and keep on ice while handling
- Do not freeze

Ordering Information

CellPrime[®] rAlbumin is provided as a 10% (w/v), 0.2 μ m filtered solution in PBS (25 mM phosphate buffer, 145 mM NaCl, pH 7.2) containing 2 and 5 g of recombinant albumin, respectively. For research and manufacturing use only.

Product Name	Volume (mL)	rAlbumin (g)	Cat. No.
CellPrime® rAlbumin	20	2	1.06331.0020
CellPrime® rAlbumin	50	5	1.06331.0050

MilliporeSigma 400 Summit Drive Burlington, MA 01803



To place an order or receive technical assistance, visit EMDMillipore.com/contactPS

For additional information, visit **EMDMillipore.com**

© 2020 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. MilliporeSigma, the vibrant M, SAFC, CellPrime and Emprove are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources. MS_PF6170EN Ver. 1.0 32060 06/2020