# Sigma-Aldrich.

Product Information

## Ribonuclease A from bovine pancreas

Sigma Type I-A, ≥60% RNase A basis (SDS-PAGE), ≥50 Kunitz units/mg protein

#### R4875

## **Product Description**

CAS Registry Number: 9001-99-4

Enzyme Commission (EC) Number: 3.1.27.5

Synonyms: RNase A, Pancreatic ribonuclease, Ribonuclease 3'-pyrimidinooligonucleotidohydrolase,

Ribonuclease I, Endoribonuclease I Molecular mass: 13.7 kDa (based on

amino acid sequence)

Extinction coefficient:  $^2$  E<sup>1%</sup> = 7.1 (280 nm)

Isoelectric point: $^{3}$  pI = 9.6

Optimal temperature: 60 °C (activity range of

15-70 °C)

Optimal pH:<sup>4</sup> 7.6 (activity range of 6-10)

Inhibitors: ribonuclease inhibitor

RNase A is an endoribonuclease that attacks at the 3'-phosphate of a pyrimidine nucleotide. For example, RNase A will cleave pG-pG-pC-pA-pG to give pG-pG-pCp and A-pG. The highest activity is exhibited with single-stranded RNA.<sup>5</sup>

RNase A is a single chain polypeptide with 4 disulfide bridges. In contrast to RNase B, RNase A is not a glycoprotein.<sup>6</sup> RNase A can be inhibited by alkylation of His<sup>12</sup> or His<sup>119</sup> (present in the active site of the enzyme).<sup>7</sup> Activators of RNase A include potassium and sodium salts.

Several theses<sup>8</sup> and dissertations<sup>9-17</sup> have cited use of product R4875 in their protocols.

### Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

## Solubility

RNase A is soluble in water at 10 mg/mL.

## **Preparation Instructions**

Solutions prepared from RNase A lyophilized powder can be made free of DNase by boiling. One procedure is as follows, according to one literature method:<sup>18</sup>

- 1. Prepare a 10 mg/mL stock solution in 10 mM sodium acetate buffer, pH 5.2.
- 2. Heat to 100 °C for 15 minutes. Allow to cool to room temperature.
- Adjust to pH 7.4 using 0.1 volume of 1 M Trizma<sup>®</sup>-HCl, pH 7.4.
- 4. Aliquot and store at -20 °C.

If RNase A is boiled at a neutral pH, precipitation will occur. When boiled at a lower pH, some precipitation may occur because of protein impurities that are present.

## Storage/Stability

This product is supplied as a lyophilized powder. Store RNase A at -20 °C.

RNase A is a very stable enzyme. Stock solutions stored in frozen aliquots remain active for at least 6 months. RNase A solutions have been reported to withstand temperatures up to 100 °C. At 100 °C, an RNase A solution is most stable between pH 2.0-4.5.<sup>19</sup>

## Usage

1

A major application for RNase A is the removal of RNA from preparations of plasmid DNA. For this application, DNase-free RNase A is used at a final concentration of 10 µg/mL.<sup>20</sup>

#### References

- 1. Smyth, D.G. et al., J. Biol. Chem., **238(1)**, 227-234 (1963).
- Keller, P.J. et al., J. Biol. Chem., 233(2), 344-349 (1958).



- 3. Tanford, C., and Hauenstein, J. D., *J. Am. Chem. Soc.*, **78(20)**, 5287-5291 (1956).
- 4. Schomberg, D., and Salzmann, M., *Enzyme Handbook*, Vol. 3, 1-3, under EC 3.1.27.5 (1990).
- Burrell, M.M., Methods Mol. Biol., 16, 263-270 (1993).
- Plummer, T.H., Jr., and Hirs, C.H.W., J. Biol. Chem., 238(4), 1396-1401 (1963).
- Heinrikson, R.L. et al., J. Biol. Chem., 240(7), 2921-2934 (1965).
- Nicolini, Anthony M., "N-Isopropyl-Acrylamide Conjugated Polyglycerol as a Delivery Vehicle for in vitro siRNA Transfection". Georgia Institute of Technology, M.S. thesis, p. 45 (2011).
- 9. Liu, Weiguo, "Transformation of Microspores for Generating Doubled Haploid Transgenic Wheat (*Triticum aestivum* L.)". Washington State University, Ph.D. dissertation, p. 96 (2004).
- Balan, Sibu, "Disulfide Bridging Poly(ethylene glycol) Reagents For Site-specific Protein Conjugation". University of London, Ph.D. dissertation, p. 71 (2007).
- 11. Michel, Dian Agustina, "Implementation and optimization of a large-scale ENU mouse mutagenesis screen and characterisation of five ENU-induced limb mutant lines". Technischen Universität München, Dr. rer. nat. dissertation, p. 55 (2009).
- 12. Tiedau, Daniela, "Tudor domain containing protein 6 and its essential role in murine spermatogenesis". Technischen Universität Dresden, Dr. rer. nat. dissertation, p. 83 (2009).
- 13. Xu, Yan, "Investigation of key non-coding and coding genes in cutaneous melanomagenesis". University of Edinburgh, Ph.D. dissertation, p. 58 (2011).
- 14. Hauff, Martha J., "An Investigation of the Nutritional Condition of Low-Latitude Fish Larvae: Growth, Transport, and Implications for Population Connectivity". University of Miami, Ph.D. dissertation, pp. 19, 50, 97 (2012).
- Minchell, Nicola E., "DNA Topological Stress during DNA Replication in Saccharomyces cerevisiae". University of Sussex, Ph.D. dissertation, p. 59 (2018).

- 16. Chibaya, Loretah, "Akt Regulation of Mdm2-p53 Signaling in Cellular Stress Responses and Tumorigenesis". University of Massachusetts Graduate School of Biomedical Sciences, Worcester, Ph.D. dissertation, p. 118 (2019).
- 17. Kumari, Poonam, "Sirt7 acts as a novel regulator of chromatin dynamics through modulation of the Sirt1/Suv39H1 axis: Possible role in cancer". Justus-Liebig-Universität Gießen, Dr. rer. nat. dissertation, p. 33 (2020).
- 18. Sambrook, J., and Russell, D.W., *Molecular Cloning, A Laboratory Manual* (3<sup>rd</sup> ed). Cold Spring Harbor Laboratory Press (Cold Spring Harbor, NY), Volume 3, A4.39 (2001).
- 19. Crestfield, A.M. et al., J. Biol. Chem., **238(2)**, 618-621 (1963).
- Sambrook and Russell, Volume 1, 1.78-1.79 (2001).

## **Notice**

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

## **Technical Assistance**

Visit the tech service page at SigmaAldrich.com/techservice.

## Standard Warranty

The applicable warranty for the products listed in this publication may be found at SigmaAldrich.com/terms.

### **Contact Information**

For the location of the office nearest you, go to <a href="SigmaAldrich.com/offices">SigmaAldrich.com/offices</a>.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

